


**CYTEC**

*Technology ahead of its time™*

[www.cytec.com](http://www.cytec.com)

# CYPHOS® IL 106

## Phosponium Ionic Liquid

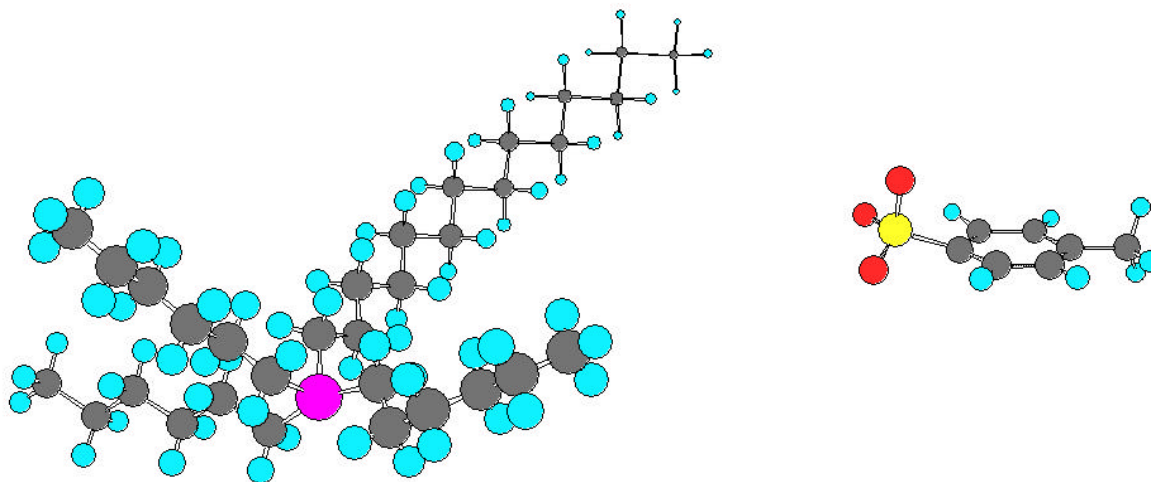

## Introduction:

### Why consider a “phosphonium” ionic liquid?

Soon after the discovery that certain nitrogen based room temperature liquid salts were found to be useful as battery electrolytes (1a,1b,), interest in these and similar salts as novel fluids and solvents developed. There were a scant number of papers during the 1980s and early 1990s but mainly due to the efforts of the group at The Queen’s University – Belfast, headed up by Professor Ken Seddon, there has been an exponential rise in interest and number of publications in the last 7 to 8 years. (2) Indeed, almost an entire issue of Green Chemistry (3) has been devoted to ionic liquids.

Perhaps one of the most influential publications to direct industrial attention to ionic liquids was a feature article entitled “Designer Solvents” in C&E News – March 30,1998 in which Ken Seddon, Robin Rogers, Tom Welton, Helene Olivier and others elaborated on the potential of ionic liquids. While the article dealt almost entirely with nitrogen based ionic liquids, there was a brief reference by Ken Seddon which alluded to the fact that phosphonium salts are also a potential source of numerous ionic liquids. This brief reference to phosphonium ionic liquids is very much representative of the current fraction of publications relating to phosphonium based ionic liquids. With the exception of several papers and patents by George Parshall in the mid 1970s using stannate and germanate salts and John Knifton et al in the early 1990s which centre on the use of molten tetrabutylphosphonium bromide as an ionic solvent, almost the entire volume of ionic liquid literature deals with nitrogen based systems and in particular, those based on 2-methylimidazolium salts.

There was a good reason for the lack of phosphonium based ionic liquid publications – availability of the starting material! While Cytec has been commercially producing phosphine derivatives since 1971, it was not until 1990 that tributylphosphine was produced on a large commercial scale. Since that time, not only has tetrabutylphosphonium chloride and bromide become available in multi ton scales, many other trialkylphosphines and the corresponding quaternary phosphonium salts are or can be manufactured on a large scale.

The phosphonium cation contains four substituents and the various combinations along with the multitude of various available anions represents an enormous number of possible salts. Even when one restricts the cation to the generic formula –  $[PR_3R']^+$ , the number is still very large. Of course, not all such phosphonium salts are liquid at room temperature, but by a judicious selection of R and R’ as well as the appropriate anion, there are many phosphonium salts which are in fact liquid at room temperature and many more which fall within the broad general definition of ionic liquids as salts which are low melting – that is less than 100 °C.

There are several reasons why one might consider a phosphonium ionic liquid. The most important one for those contemplating an industrial process is availability and cost. Phosphonium salts can meet both of these demands – already Cytec is manufacturing phosphonium salts on a multi ton scale and because of the high volumes, costs will be relatively low. For commercial products, chemical inventory registration is also part of the availability equation. While, most of the possible phosphonium ionic

liquids are still not registered, several are already listed on EINECS, TSCA, EEC, AICS, PICCS and DSL.

Ionic liquids, in general, are not going to be outrageously expensive, but they will not be in the same league as toluene, 2-hydroxypropane ( IPA ) or tetrahydrofuran ( THF ). This means that to be economically viable, they must be chemically as well as thermally very stable for multiple recycle use. Even 0.5 to 1% decomposition can lead to major losses after 10 to 20 cycles. Not only will there be solvent losses but there will also be contamination of the ionic liquid solvent and/or products with decomposition byproducts.

In this regard, phosphonium salts are much more thermally stable than the corresponding ammonium salts and even have an edge on imidazolium salts. This is very important for processes which operate at temperatures greater than 100 °C. In addition to being slightly less thermally stable, the imidazolium cation contains protons which are not entirely inert. They are some what acidic which can result in carbene formation. Phosphonium salts, on the other hand have no such acidic protons.


The fact that alkylphosphonium salts are, in general, less dense than water can be beneficial in product work-up steps while decanting aqueous streams which contain inorganic salt byproducts. Imidazolium salts, on the other hand are more dense than water.

**Trade Name :** **CYPHOS IL 106 phosphonium ionic liquid**

**Chemical Name:** triisobutyl(methyl)phosphonium tosylate

**C.A.S. Number:** None

**Registration:** None

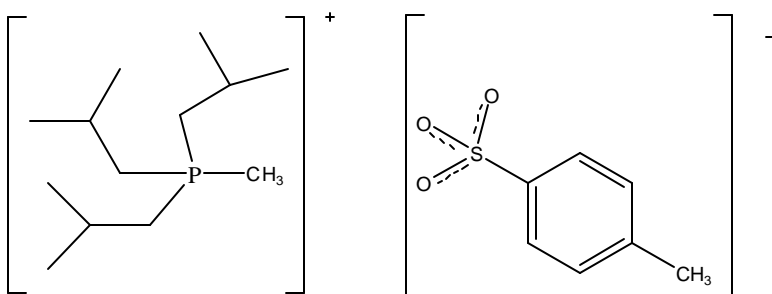

**CYPHOS IL 106 phosphonium ionic liquid** is prepared from triisobutylphosphine and methyltosylate and consequently is **totally halogen free**. ( 3,8 )

## Density:

A plot of density as a function of temperature for **CYPHOS IL 106** can be found in Figure 1. The presence of the aromatic anion results in an increased density and consequently, **CYPHOS IL 106** is one of the few phosphonium ionic liquids with a density greater than that of water.

## Miscibility:

**CYPHOS IL 106** is totally miscible with water as well as with common organic solvents such as those listed in Table 1.

**Table 1**

### **CYPHOS IL 106 Miscibility**

<b>Diluent</b>	<b>Miscible</b>
water	Yes
hexane	No
toluene	Yes
2-hydroxypropane (IPA)	Yes
diethylether	Yes
tetrahydrofuran	Yes
dichloromethane	Yes

## Viscosity:

It is a moderately viscous liquid with a density of approximately 1. ( Figure 2 ). However, as with all ionic liquids, even the addition of the slightest amount of solute will have a dramatic effect on lowering the viscosity. ( 8 ), ( Figure 3 ). The viscosity also decreases exponentially with temperature ( Figures 2 and 3 ), so at typical reaction temperatures of 80 to 100 °C, the **CYPHOS IL 106** system will become quite water-like in viscosity.

## Thermal Stability:

Typically standard TGA plots are used to determine relative thermal stability of ionic liquids. Heating rates of 5 to 10 °C per minute either under an inert atmosphere or under oxidative conditions such as air are usually reported. Under these conditions, the onset for weight loss for **CYPHOS IL 106** is approximately 320 and 380 °C under dinitrogen and air respectively. ( Figure 4 ). However, in reality, the true temperature at which an ionic liquid is thermally stable is much lower. Figures 5 and 6 are isothermal TGA plots under dinitrogen and air respectively. The respective safe operating temperatures would appear to be 180 and 210 °C.

## Applications:

Phosphonium tosylates have been reported as novel solvents for the hydroformylation of olefins. ( 7 )

Bio-transformations are often limited to the concentration of either the substrate or products due to toxicity effects on either the enzymes or organisms. For example, 1500 mg/l of phenol is toxic to *Pseudomonas putida*. However, the bio-degradation can be carried out in a two phase system such as **CYPHOS IL 106** / (aqueous phenol), in which the phenol will partition into the IL phase at a concentration which is non-toxic to *P. putida*. This organism is totally compatible with **CYPHOS IL 106** and is thus able to consume the phenol from the aqueous phase as it gradually partitions into the IL phase (9).

## Analysis:

An assay procedure for **CYPHOS IL 106** is currently under development. However, based on  $^{31}\text{P}$  NMR data, it is expected to assay > 95%. The main impurities are likely to be triisobutylphosphine oxide and or a trace of methyltosylate. Water may also be present to the extent of 0.2%, but the important feature is that **CYPHOS IL 106** will be totally halogen free.

$^{31}\text{P}$  NMR is generally not suitable for precise analysis. However, the distinctive signal at +33 ppm can be useful for qualitative analysis.

While electrospray mass spectral analysis ( ESMS ) is generally not available in every laboratory, this also a very useful tool for both quantitative and qualitative analysis.

LC/MS is the most universal tool to identify and quantify cations and anions (10).

## References:

- 1a) Fannin, A .A, Jr; King, L. A.; Levisky, J. A.; Wilkes, J. S.; J. Phys. Chem., 1984, **88**, 2610-2614
- 1b) Fannin, A. A. Jr; Floreani, D. A.; King, L. A.; Landers, J. S.; Piersma, B. J.; Stech, D. J.; Vaughn, R. L.; Wilkes, J. S.; and Williams, J. L., J. Phys. Chem. 1984, **88**, 2614-2621
- 2) "Forward", Professor Ken Seddon, Green Chemistry April, 2002.
- 3) Green Chemistry April, 2002

- 4) "Industrial Preparation of Phosphonium Ionic Liquids"; Christine Bradaric, Andrew Downard, Christine Kennedy, Allan Robertson, Yuehui Zhou; Green Chemistry, 2003, **5**, 143-152
- 5) "Phosphonium Salts"; A Robertson; WO 01/87900 A1
- 6) "Process for Preparing Ambient Temperature Ionic Liquids"; Kenneth Richard Seddon, Adrian John Carmichael, Martyn John Earle; WO0140146
- 7) "Clean catalysis with ionic solvents-phosphonium tosylates for hydroformylation"; Karodia, Nazira; Guise, Steven; Newlands, Craig; Andersen, Jo-Ann; Chemical Communications (Cambridge) (1998), (21), 2341-2342
- 8) K.R. Seddon, A. Stark, M.J. Torres; Pure Appl. Chem. , 2000, **72**, 2275
- 9) "Phosphonium ionic liquids for degradation of phenol in a two-phase partitioning bioreactor"; Environmental Biotechnology; 2005, *67*, 131-137; M.D. Baumann, A.J. Daugulis and P.G. Jessop
- 10) "Development of an LC/MS Analytical Procedure for Ionic Liquids with Emphasis on Phosphonium Ionic Liquids"; 1<sup>st</sup> International Congress on Ionic Liquids, June 19-22, 2005, Salzburg, Austria; T.T. Chang and M.J. Piquette ( also posted on [www.cyttec.com/business/phosphine](http://www.cyttec.com/business/phosphine) )

Figure 1

**CYPHOS IL 106 – Density vs. Temperature**

CYPHOS IL 106 - Density vs. Temperature

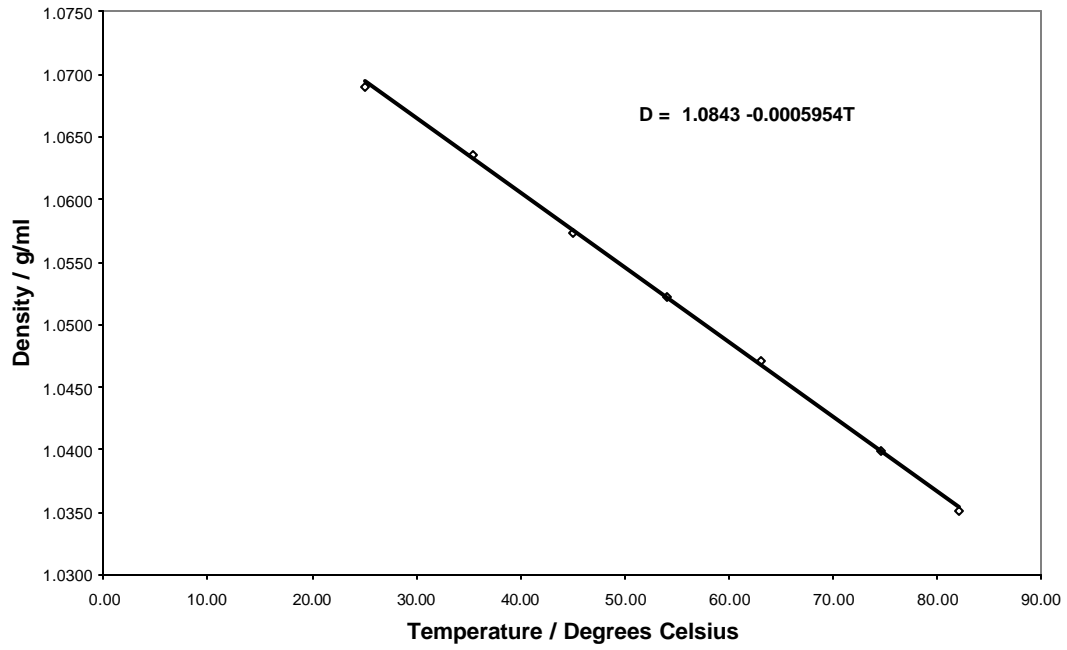


Figure 2

**CYPHOS IL 106 – Viscosity vs. Temperature**

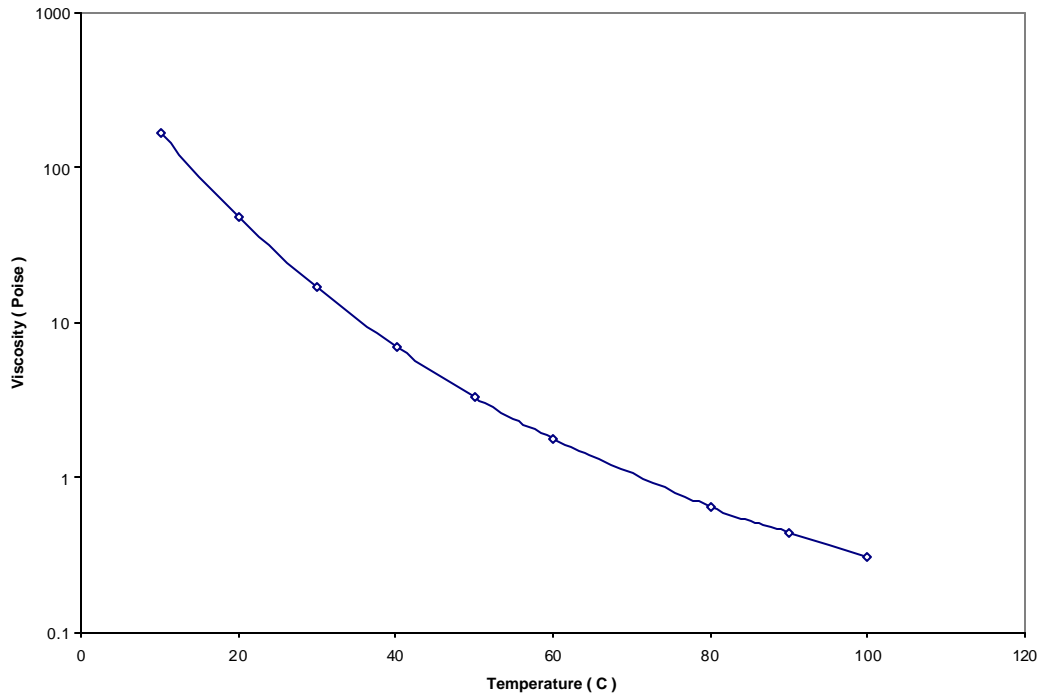


Figure 3

**CYPHOS IL 106 – Effect of Water and Toluene on Viscosity**


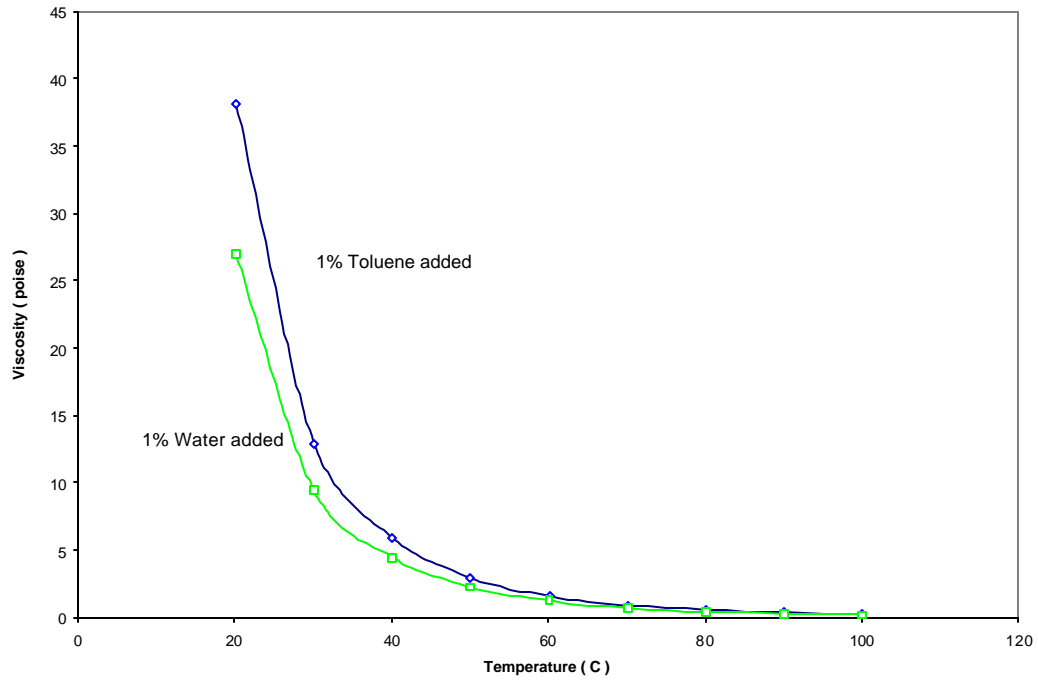



Figure 4

**CYPHOS IL 106 – Standard TGA Plots under Air and N<sub>2</sub>**

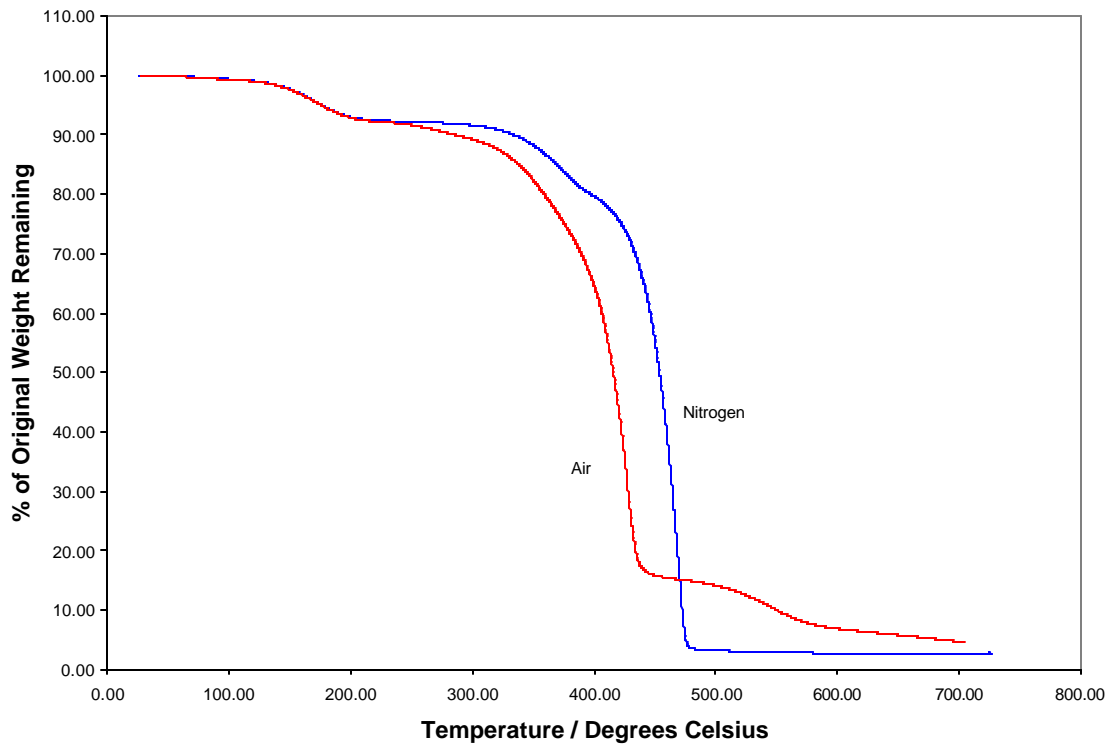


**Figure 5**

**CYPHOS IL 106 – Isothermal TGA Plots under Air**

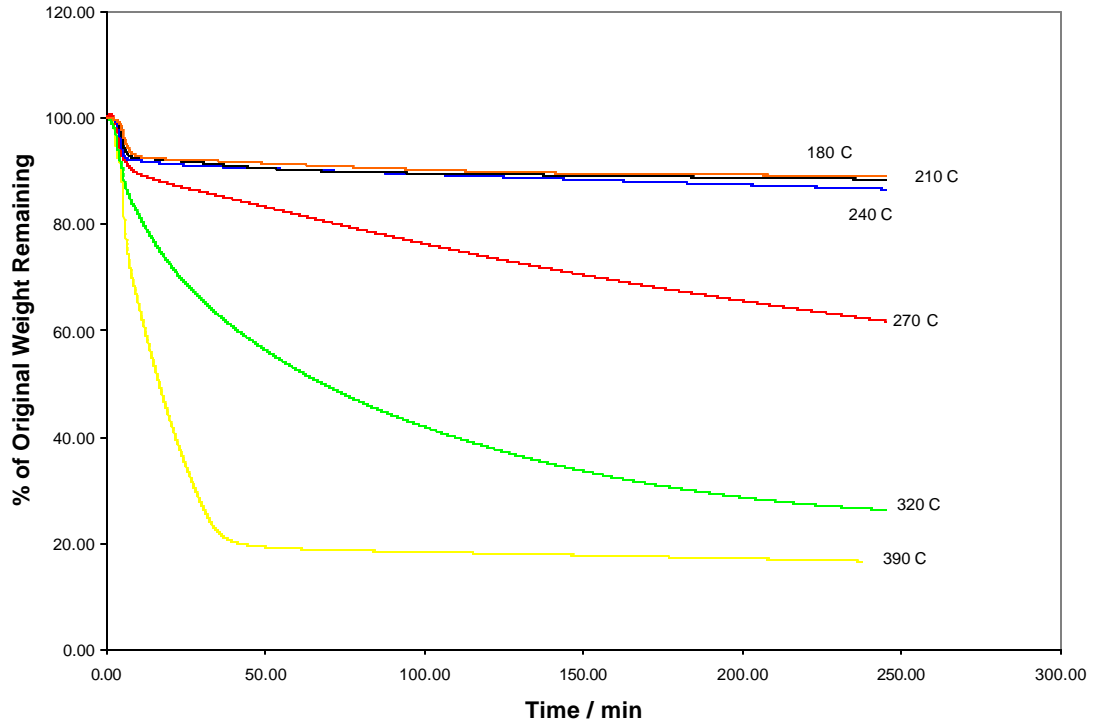

**Figure 6**

**CYPHOS IL 106 – Isothermal TGA Plots under N<sub>2</sub>**

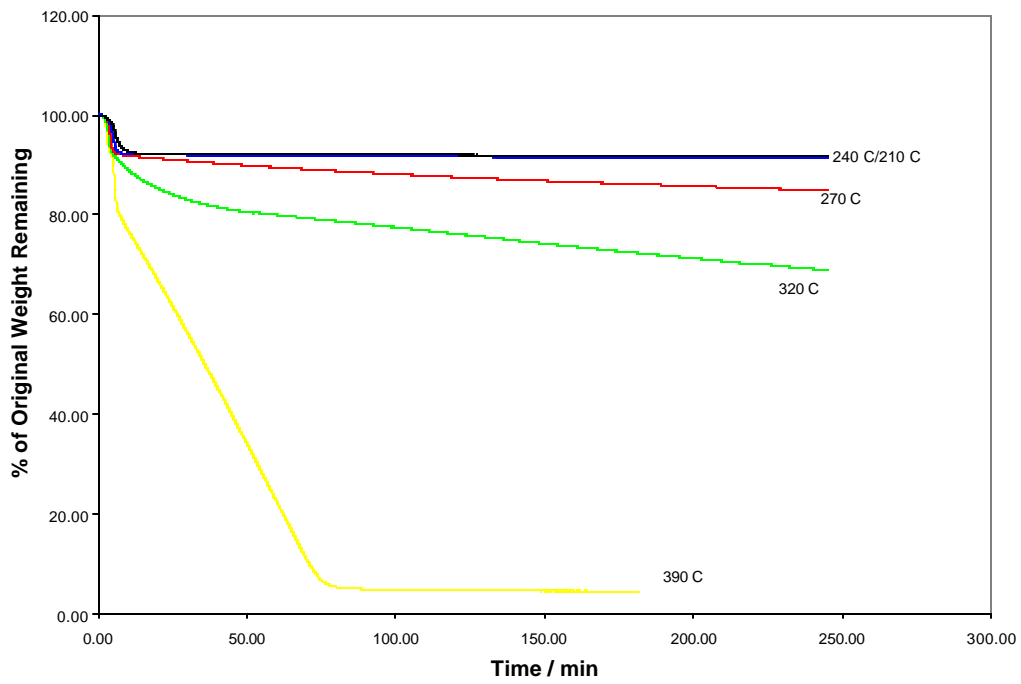

**Important Notice:** The information and statements in this data sheet are believed to be reliable but are not to be construed as a warranty or representation for which we assume legal responsibility or as an assumption of a duty on our part. Users should undertake sufficient verification and testing to determine the suitability for their own particular purpose of and information or products or vendors referred to within. **NO WARRANTY OF FITNESS FOR A PARTICULAR PURPOSE IS MADE.** Nothing in this data sheet is to be taken as permission, inducement or recommendation to practice any patented invention without a license.

**Trademark Notice:** The ® indicates a Registered Trademark in the United States and the TM or \* indicates a Trademark in the United States. The mark may also be registered, the subject of an application for registration or a trademark in other countries.

**Last updated: September 2, 2005**